



Check veterinary drug & pesticides in honey with **Ringbio Rapid Test Kit**

- 🌸 One Step Universal Preparation
- 🌸 Rapid 10min Test
- 🌸 No Instrument Required



RING BIOTECHNOLOGY CO LTD



Scan to read PDF copy

Drug Residue Rapid Test Kit

Item #	Product - 48 tests/kit
100307	Tetracyclines Rapid Test Kit
100309	Quinolones Rapid Test Kit Kit
100314	Chloramphenicol Rapid Test Kit
100323	Tilmicosin & Tylosin, Macrolides Rapid Test Kit
100334	Carbendazim Rapid Test Kit
100335	Acetamiprid Rapid Test Kit
100336	Spectinomycin, Thiamphenicol & Quinolones, SQT Rapid Test Kit
100337	Tetracyclines, Macrolides & Erythromycin, TME Rapid Test Kit
100338	Carbendazim, Acetamiprid & Carbaryl, MDL Rapid Test Kit
100339	Imidacloprid, Cyromazine, Thiamethoxam, Carbofuran, DEMN Rapid Test Kit

We also have ELISA kit for honey testing. Contact for details.

Procedures for testing



Take honey into a tube and mix with sample buffer



Add 200ul mixed sample into vial mix and incubate at 40°C Wait for 5min



Insert test strip Wait for 5min at 40°C



Check the result

** this procedure applies for drug residue with MRL.

For not allowed drugs such as nitrofurans, please contact us



Ring Biotechnology Co Ltd

Add: Building 3, Zhongtongtai TechnoPark, BDA, Beijing 100176, China

Tel: +86-10-56267496 / +86-10-26242030

Email: info@ringbio.com

Web: www.ringbio.com