




Check the meat and shrimp with Antibiotic Residue Rapid Test Kit

- 🌻 One Step
- 🌻 Rapid 10min Test
- 🌻 No Instrument Required



RING BIOTECHNOLOGY CO LTD

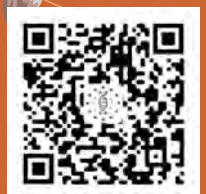


Scan to read PDF copy

Antibiotic Residue Rapid Test Kit

| Item # | Product - 48 tests/kit | Sample |
|--------|--|--------|
| 510001 | Meat Beta Lactams Rapid Test Kit | Meat |
| 510005 | Meat Lincomycin Rapid Test Kit | Meat |
| 510006 | Meat Tetracyclines Rapid Test Kit | Meat |
| 510008 | Meat Sulfonamides Rapid Test Kit | Meat |
| 510009 | Meat Fluoroquinolones Rapid Test Kit | Meat |
| 510014 | Meat Chloramphenicol Rapid Test Kit | Meat |
| 510023 | Meat Macrolides Rapid Test Kit | Meat |
| 510032 | Meat Florfenicol Rapid Test Kit | Meat |
| 510034 | Meat Trimethoprim Rapid Test Kit | Meat |
| 510035 | Meat Olaquinox Rapid Test Kit | Meat |
| E4701 | Nitrofurantoin AOZ/AMOZ/AHD/SEM 4-in-1 QuaTest | Shrimp |
| E1705 | Chloramphenicol Rapid Test Kit | Shrimp |

We also have ELISA Kits for meat and shrimp testing. Contact us for details.


Procedures for testing

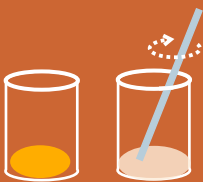

Homogenized meat into a tube and mix it with sample buffer



Centrifuge

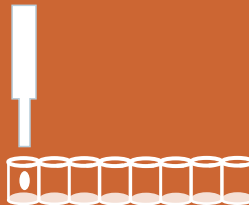

Add 200ul transparent layer into vial mix and wait for 5min



Insert test strip
Wait for 5min
Check the result

** this procedure applies for drug residue with MRL.

For not allowed drugs such as CAP and nitrofurantoin, please contact us



Ring Biotechnology Co Ltd

Add: Building 3, Zhongtongtai TechnoPark, BDA, Beijing 100176, China

Tel: +86-10-56267496 / +86-10-26242030

Email: info@ringbio.com

Web: www.ringbio.com